

## Operating Instructions




### OIV 255 Software-Update to V3.16.0.0

#### Preparing the data

At first, the USB stick should be insert into the OIV-2 so that the microscope system creates the corresponding folder. Then remove the stick and connect it to a PC. After that, copy the zipped file (*UpdateMeasure\_R\_SYGD\_V3.16.0.0\_20210531.tar.gz*) and past it in the **update** folder (Path: USB disk/camera/update). Finally, insert the USB stick into OIV-2 again, and upgrade according to the previous operation method.


#### Operating method

Auxiliary tools > Advanced setting


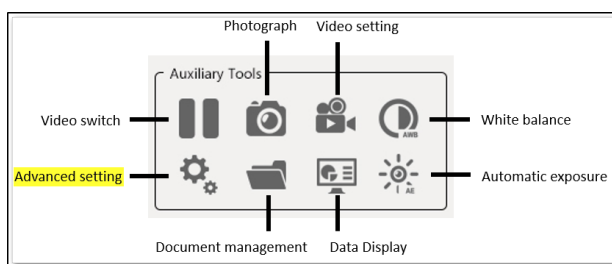


## Operating Instructions

### Setting > Update

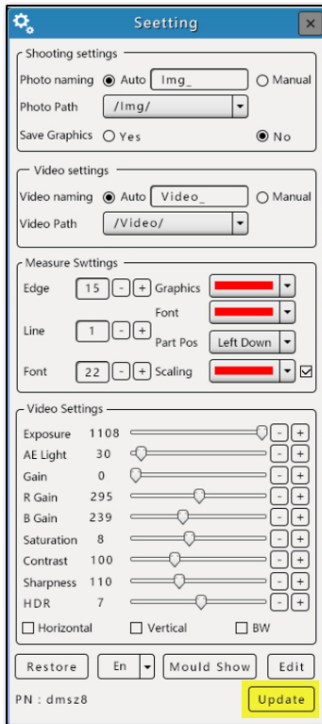

Select the file (Package name) > Update > OK

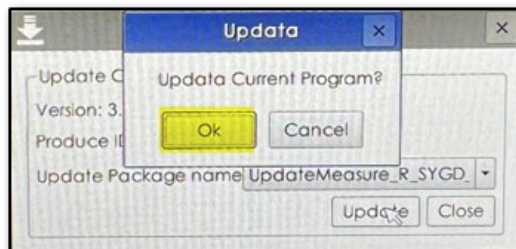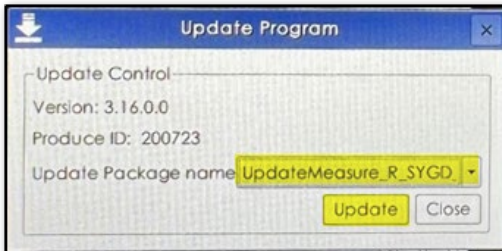

After the update has been finished, the language must be adapted too. For that go back to Auxiliary tools > Setting and select EN

